

TECHNICAL DATA SHEET

Article No. 9721

Tryptic Soy Broth (TSB) Ph. Eur.

SYNONYMS

CASO Bouillon, Casein Soya bean Digest Broth

SPECIFICATION

Highly nutrient liquid medium for general purpose use, formulated according to Pharmacopeial Harmonized Method.

FORMULA* IN G/L

Casein peptone	17.0
Soy peptone	3.0
Sodium chloride	5.0
Dipotassium phosphate	2.5
Dextrose	2.5

Final pH 7.3 ±0.2 at 25 °C

*Adjusted and/or supplemented as required to meet performance criteria.

DIRECTIONS

Dissolve 30 g of powder in 1 l of distilled water. Distribute in suitable containers and sterilize by autoclaving at 121 °C for 15 minutes.

DESCRIPTION

The Tryptic Soy Broth was initially developed for the cultivation of very fastidious microorganisms without the addition of serum, blood or any other enrichment agent.

As a general purpose culture medium it supports the growth of most organisms, both aerob and facultative anaerobes, even if their requirements are high. Due to its high vitamin content *Brucella*, *Pasteurella* and *Streptococcus* are perfectly viable, moreover a CO₂ enriched atmosphere can further enhance growth.

In anaerobic conditions this broth will grow *Bacteroides* and *Clostridium* species. For this purpose, the best results can be obtained by adding 0.3 % agar and 0.05 % sodium azide for *Clostridium*.

Th. Geyer GmbH & Co. KG

Dornierstr. 4 – 6
D-71272 Renningen
Tel.: +49 7159 1637-0
Fax: +49 7159 1637-710
renningen@thgeyer.de
www.thgeyer.de

BW-Bank (Swift/BIC SOLADEST600)
IBAN DE85600501010002036302
Postbank Stuttgart (Swift/BIC PBNKDEFFXXX)
IBAN DE3260010070000020708
Deutsche Bank (Swift/BIC DEUTDESSXXX)
IBAN DE06600700700125518100

St.-Nr. 70093/40018 / USt-IdNr. DE147510304
Amtsgericht Stuttgart / HRA-Nr. 254140
Persönlich haftende Gesellschafterin:
Geyer Beteiligungsgesellschaft mbH
Amtsgericht Stuttgart / HRB-Nr. 252035
Geschäftsführer: Lutz-Alexander Geyer / Thomas Roth

Tryptic Soy Broth's superior growth-promoting properties make it particularly suitable for tube dilution methods for antibiotic sensitivity testing.

The broth can be used for bile solubility testing in pneumococci, and also used for catalase and coagulase assays and for the preparation of hypersaline broths.

It is a most suitable medium for the preparation of antigens and toxins in bacteria, moulds and yeasts.

TSB is used as a primary enrichment medium for food examination. In the dairy industry it is employed for testing resazurine reduction.

The medium is not suitable for maintenance purposes since carbohydrate fermentation liberates many acids which may threaten the organism's viability. Therefore, though it allows the growth of streptococci and *Neisseria*, these species tend to die if repeatedly sub-cultured in this medium. Such fastidious organisms are best maintained on Cystine Tryptone Fluid Medium (CTA) or even TSA (Art. no. 9738).

QUALITY CONTROL

- Incubation temperature: 30-35 °C
- Incubation time: 24-72 h
- Inoculum: Practical range 10-100 CFU. Min. 50 CFU (productivity), according to Ph. Eur.

Microorganism	Growth	Remarks
<i>Bacillus subtilis</i> ATCC® 6633	Good	None
<i>Staphylococcus aureus</i> ATCC® 6538	Good	None
<i>Escherichia coli</i> ATCC® 8739	Good	None
<i>Pseudomonas aeruginosa</i> ATCC® 9027	Good	None
<i>Salmonella typhimurium</i> ATCC® 14028	Good	None
<i>Aspergillus brasiliensis</i> ATCC® 16404	Good	≤5 d
<i>Candida albicans</i> ATCC® 10231	Good	≤5 d

REFERENCES

- ATLAS, R.M. & L.C. PARKS (1993) Handbook of Microbiological Media. CRC Press, Inc. London.
- DOWNES, F.P. & K. ITO (2001) Compendium of Methods for the Microbiological Examination of Food, 4th ed. ASM. Washington. DC.
- EUROPEAN PHARMACOPOEIA 8.0 (2014) 8th ed. § 2.6.13. Microbiological examination of non-sterile products: Test for specified microorganisms. Harmonised Method. EDQM. Council of Europe. Strasbourg.
- FDA (Food and Drug Administrations) (1998) Bacteriological Analytical Manual. 8th ed. Revision A. AOAC International. Gaithersburg. MD.
- HORWITZ, W. (2000) Official Methods of Analysis of AOAC INTERNATIONAL. 17th ed. Gaithersburg. MD. USA.

Th. Geyer GmbH & Co. KG

Dornierstr. 4 – 6
D-71272 Renningen
Tel.: +49 7159 1637-0
Fax: +49 7159 1637-710
renningen@thgeyer.de
www.thgeyer.de

BW-Bank (Swift/BIC SOLADEST600)
IBAN DE85600501010002036302
Postbank Stuttgart (Swift/BIC PBKDEFFXXX)
IBAN DE32600100700000020708
Deutsche Bank (Swift/BIC DEUTDESSXXX)
IBAN DE06600700700125518100

St.-Nr. 70093/40018 / USt-IdNr. DE147510304
Amtsgericht Stuttgart / HRA-Nr. 254140
Persönlich haftende Gesellschafterin:
Geyer Beteiligungsgesellschaft mbH
Amtsgericht Stuttgart / HRB-Nr. 252035
Geschäftsführer: Lutz-Alexander Geyer / Thomas Roth

- ISO 11133:2014. Microbiology of food, animal feed and water. Preparation, production, storage and performance testing of culture media.
- ISO 9308-1 Standard (2000) Water Quality. Detection and enumeration of E. coli and coliform bacteria. Membrane filtration method.
- PASCUAL ANDERSON, M^ªR^a (1992) Microbiología Alimentaria. Díaz de Santos S.A., Madrid.
- USP 33 - NF 28 (2011) <62> Microbiological examination of non-sterile products: Test for specified microorganisms. Harmonised Method. USP Corp. Inc. Rockville. MD. USA.
- USP 33 - NF 28 (2011) <71> Sterility Tests. Harmonised Method. USP Corp. Inc. Rockville. MD. USA.

STORAGE

Keep tightly closed, away from light, in a dry place (4-30 °C).

SHELF LIFE

5 years from date of production.

Th. Geyer GmbH & Co. KG

Dornierstr. 4 – 6
D-71272 Renningen
Tel.: +49 7159 1637-0
Fax: +49 7159 1637-710
renningen@thgeyer.de
www.thgeyer.de

BW-Bank (Swift/BIC SOLADEST600)
IBAN DE85600501010002036302
Postbank Stuttgart (Swift/BIC PBNKDEFFXXX)
IBAN DE32600100700000020708
Deutsche Bank (Swift/BIC DEUTDESSXXX)
IBAN DE06600700700125518100

St.-Nr. 70093/40018 / USt-IdNr. DE147510304
Amtsgericht Stuttgart / HRA-Nr. 254140
Persönlich haftende Gesellschafterin:
Geyer Beteiligungsgesellschaft mbH
Amtsgericht Stuttgart / HRB-Nr. 252035
Geschäftsführer: Lutz-Alexander Geyer / Thomas Roth